



— Stock Code: 688026 —



— Stock Code: 688026 —

YOUR RELIABLE PARTNER IN LIFE SCIENCE



Guangzhou Jet Bio-Filtration Co.,Ltd.

Address: No.1 DouTang Road, YongHe Development Zone, Guangzhou, 511356,China
Contact Hotline: +86-400-8717-688
Before-sale service: info@jetbiofil.com
After-sale service: info@jetbiofil.com

• JET · Europe Office

Akralab, S.L.
Pol. Ind. Las Atalayas, Av. de la Antigua Peseta, 77,
Buzón 20212, 03114 Alicante, Spain Buzón 20212
Tel:+902 22 22 75 | +34 965 11 65 21
E-mail:mariano.guzman@akralab.es www.akralab.es

• JET · Israel Office

18 Soreq St. Emek Sara 8400795,
Beer-sheva, Israel
Tel: 972-86280210 Fax:972-8-6280801
E-mail: info@romical.com

• JET · Japan Office

Manshon Omiya 206, Mibuai-cho 32,
Nakagyo-ku, Kyoto 604-8812, Japan
Tel:+81-75-334-6282 Fax:+81-75-822-2194
E-mail:kooki@ic5.so-net.ne.jp

• JET · Korea Office

DAIHAN Bldg.,24-4, Sangwolgok-dong,Seongbuk-gu,
Seoul,Korea
Tel: +82-80-008-3000 Fax:+82-33-737-7509
E-mail: chloe.baek@daihan-sci.com

• JET · South America Office

Gral. Alvear 123, 1° Of. 5 Lomas de Zamora (B1832BVC),
Buenos Aires, Argentina
Tel: (+5411) 42921611
E-mail: info@apbiotech.com.ar www.apbiotech.com.ar

• JET · Australia Office

11 Challenge Boulevard, Wangara WA 6065 Australia
Tel: +61 8 9302 1911
E-mail: rowewa@rowe.com.au

• JET · US Office

1551 South Acottdale Court Suirt 200,Elgin IL,60123
Tel: 847-622-0456
E-mail: pkne007@aol.com



Pipette Micro Tips Technical Manual



Guangzhou Jet Bio-Filtration Co.,Ltd.





Contents

Abstract

Molecular test is a laboratory test that is used to study constituent cells and body fluids using DNA and/or RNA detection technology to identify the molecular characteristics and abnormalities under the basic principle of PCR. Molecular tests are widely used in various fields, such as laboratories, and clinical and non-clinical fields. Molecular diagnosis, an example of the application of molecular tests for in vitro diagnosis, has currently become the fastest-growing and cutting-edge technology in the field of in vitro diagnosis. In addition to disease diagnosis, scientific research institutes, pharmaceutical companies, and CROs also use molecular test technologies and products to carry out research and development.

Consumables for molecular tests produced by Jet Biofil are DNase/RNase and pyrogen-free and produced in Class 100,000 cleanrooms with high-quality raw materials conforming to USP Class VI standards. The pipette tips have a variety of specifications, allowing them to be compatible with various automatic instruments.

• Company Introduction	2
• Qualify Certification & Benchmarks	5
• Facilities	6
• Pipette Product Series	9
Pipette Micro Tips	9
ZEROTIP® Pipette Micro Tips	14
• Product Datasheet	18
Pipette Micro Tips	18
ZEROTIP® Pipette Micro Tips	23
• Raw Material Certifications	28
• Product Certifications of Quality	33
Pipette Micro Tips	33
ZEROTIP® Pipette Micro Tips	34
• After-Sale Supports	35
• Pipette Tips Adapter Compatibility List	36